## **CHLORINE SANITISING TABLETS**

## 3.2g Effervescent Chlorine Tablet Dilution Guidelines

DISINFECTION TASK	AVAILABLE Cl. (ppm)	WATER FOR 1 TABLET	COMMENTS
Blood spillages	10,000	100ml	Always use gloves. Cover spill liberally. Leave 2 mins before safely removing and disposing
Concrete porous floor	5,000	200ml	Scrub thoroughly with solution, remove excess and allow to air dry
Lab discard jars etc.	2,500	400ml	Immerse overnight
Environmental disinfection	1,000	1L	Immerse overnight
General lab disinfection	1,000	1L	Saturate clean cloth, wipe surfaces and equipment. Leave 5 mins min and allow to air dry or use paper towels
Stainless steel instruments	600	1.6L	Clean first then soak for 1 hour. Do not rinse.
WCs, drains	400	2.5L	Pour in at night or during quiet periods
Food prep surfaces	200	5L	After cleaning, use solution to disinfect. Leave 3 mins, air dry or use disposable paper towels
Baby bottles, teats etc.	125	8L	Thoroughly clean first. Soak for 30 mins. Rinse before use with pre boiled water
Dishcloths and Mops	60	16L	After rinsing, soak for 30-60 mins. Do not soak overnight
Salad	100	10L	Soak for 10 mins, rinse gently
Swimming Pools	1.0—3.0	500L	Dissolve 1-3 tablets in 1L tap water. Pour in at night or quiet periods by spreading.
Saunas & Jacuzzis	2.0—4.0	500L	Dissolve 1-3 tablets in 1L tap water. Pour in at night or quiet periods by spreading.





## **PACKAGING**

Order Code : XCST

Pack Size: 6 x 200 tablets



## **BIO PRODUCTIONS LTD**

72 Victoria Road, Burgess Hill, W Sussex, RH15 9LH T: 0044 1444 244000 | F: 0044 1444 244999 E: sales@bio-productions.com

W: www.bio-productions.com